

INNOVATOR INSIGHT

Scalable upstream process development for the suspension-based production of lentiviral vectors for CAR T cell therapies with multiparallel & benchtop bioreactor systems & DoE methodology

Diana Riethmüller, Alengo Nyamay'antu & Franziska Bollmann

Cell and gene-based therapies present a new treatment paradigm that have the potential to address unmet clinical needs. Viral vectors such as adenoviruses, adeno-associated viruses, retroviruses and lentiviruses are effective delivery systems for genetic material used in cell and gene therapies and vaccines. HEK-293 cells and derivatives are commonly used as a workhorse cell line for lentiviral vector (LVV) production for cell and gene therapy applications. Adherent production processes with these cells utilize static flask cultures, and this adherent method is quite easy to develop and perform. However, it also significantly lacks the ability for automation and scalability. Typical bioreactors based either on a rocking motion or stirred tank agitation can provide these features. Therefore, to scale up viral vector production for commercialization, adherent processes should be shifted to a suspension-based process, a significant challenge for the regenerative medicine industry. The use of a suspension adapted HEK-293 cell line and the Ambr[®] 15 microbioreactor system can facilitate transition from adherent cultures to suspension cultures by enabling fast process optimization with the ability to screen in parallel many parameters in small volumes. As a proof-of-concept study, we established here such a transitional protocol for the cultivation

of suspension adapted HEK-293T cells and the production of CD19-CAR lentivirus in small and benchtop scale stirred bioreactors.

Cell & Gene Therapy Insights 2021; 7(6), 689–700

DOI: 10.18609/cgti.2021.099

Cell and gene-based therapies present a new treatment paradigm that have the potential to address clinical needs that are unmet by current small molecule and biotherapeutic approaches [1].

Viral vectors such as adenoviruses, adeno-associated viruses and retroviruses are effective delivery systems for genetic material used in cell and gene therapies and vaccines. Lentiviruses are used for example for the transfer of genetic information for novel cellular immunotherapies (gene modified cell therapies), like CAR-T cell therapy [2]. These innovative approaches will be a substantial part of next-generation therapies to cure devastating diseases.

The number of clinical candidates is growing rapidly and commercial scale manufacturing is becoming a reality for these clinical candidates. The processes to manufacture viral vectors for commercialization require a high level of operator expertise and GMP guideline application; yet they are currently mainly based on an R&D approach.

HEK-293 cells and derivatives are commonly used as a workhorse cell line for lentiviral vector (LVV) production for cell and gene therapy applications. Adherent production processes with these cells utilize static flask cultures, like T-flasks, cell factories or cell stacks. This adherent method is quite easy to develop and perform. However, it also significantly lacks the ability for automation and scalability. Typical bioreactors based either on a rocking motion or stirred tank agitation can provide these features. Therefore, to scale up viral vector production for commercialization, adherent processes should be shifted to a suspension-based process, a significant challenge for the regenerative medicine industry

[3]. Suspension-based lentivirus production could either be performed using microcarriers for culturing adherent cell lines or by using a suspension adapted cell line [4]. The use of a suspension adapted HEK-293 cell line and the Ambr® 15 microbioreactor system can facilitate transition from adherent cultures to suspension cultures by enabling fast process optimization with the ability to screen in parallel many parameters in small volumes [5]. As a proof-of-concept study, we established here such a transitional protocol for the cultivation of suspension adapted HEK-293T cells and the production of CD19-CAR lentivirus in small and benchtop scale stirred bioreactors.

BENCHTOP BIOREACTOR SYSTEMS FOR SEAMLESS SCALE-UP

Ambr® 15 from Sartorius is an automated micro-scale bioreactor system that enables a fast screening of process parameters such as pH, DO, temperature, stirring rate in less time, with reduced reagents/media use and labor costs. This is a financial benefit for CDMOs and start-up companies. Parallel processing, automation capability and consistency provided by Ambr® 15 system [6,7] enables rapid, high throughput process improvement and optimization, including Design of Experiment (DoE) studies. It releases efforts for time spent in data analysis, thanks to its integration to the DoE software MODDE®.

The “big brother” of Ambr® 15, the Ambr® 250 Modular, is a bioreactor system with working volumes going from 100–250 mL and a system for scale-down model for larger stirred bioreactor systems [8]. Similarly, to the

Ambr® 15, the Ambr® 250 Modular facilitates upstream process development with reduced effort due to its parallel cultivation capacity and the possibility for “hands-off” workflow automation. The Ambr® 15 and Ambr® 250 have been shown to be valuable scale-down model systems [8,9]. Although some effort is needed to characterize them and also novel scale-down model criteria might need to be established (reviewed by [10]), due to the high throughput screening capabilities of both instruments, this characterization can be efficiently performed. Some process characteristics cannot be mimicked with the Ambr® 15 and the Ambr® 250 Modular like any continuous manipulation, e.g. perfusion or feeding cultivation strategies. Figure 1 highlights the beforementioned capabilities of the Ambr® bioreactor systems. Due to their scalability to larger stirred bioreactor systems [11], they are the ideal tools for process development.

and percentage of DO have been optimized prior to the study presented here. Secondly, we optimized the transient transfection protocol with PEIpro® transfection reagent to reach high and robust recombinant lentivirus titers. Optimization of transient transfection requires taking into account several parameters, including selection of synthetic culture media, DNA amount, ratio of DNA / PEIpro® transfection reagent and ratio of plasmids. Thirdly, we wanted to assess scalability of optimal conditions by producing lentivirus in larger volumes in the Ambr® 250 Modular. The methods used for lentivirus production and quantification were mainly based on those described by Labisch *et al.* [12]. To get meaningful results with reduced sample number, we performed a DoE study to identify optimal culture and transfection conditions by using the MODDE® software for experimental planning and analysis of results.

EFFICIENT OPTIMIZATION STRATEGY FOR SCALABLE LENTIVIRUS PRODUCTION

We first focused on optimization of cell culture conditions. We used the Ambr® 15 microbioreactor system to screen for highest viable cell count and highest lentivirus titer while varying key parameters such as stirring speed and pH value. The factors have mainly been selected based on previous findings. Some factors, e.g. the seeding cell density

OPTIMIZATION OF CULTURE CONDITIONS FOR LENTIVIRUS PRODUCTION IN AMBR® 15

To quickly identify optimal culture conditions to produce lentiviral vector, parallel analysis in an automated fashion was performed in the small-scale bioreactor Ambr® 15 system combined with DoE principles. We optimized the stirring speed and the cell culture pH within a defined range, as shown in Table 1.

► **FIGURE 1**

Sartorius stirred bioreactor portfolio for different stages of biotherapeutic manufacturing.

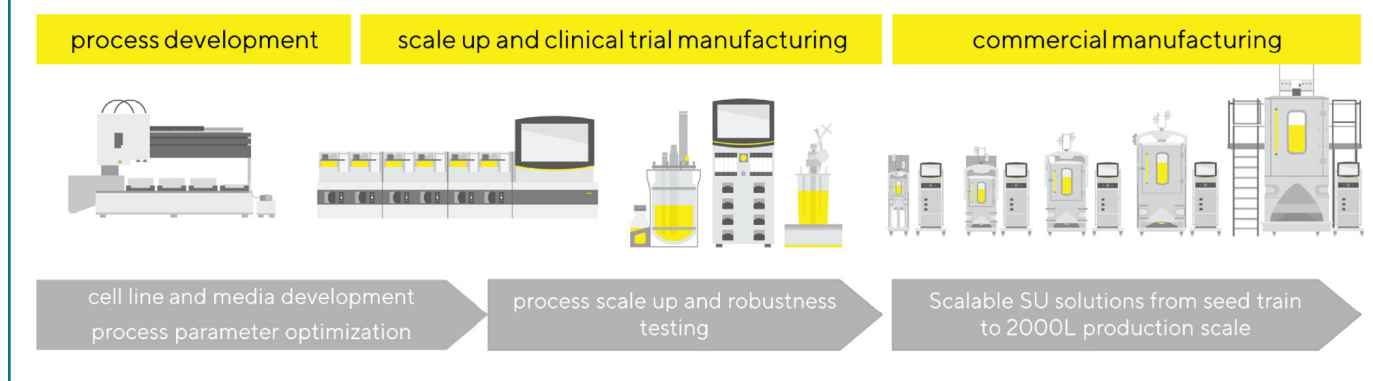


TABLE 1
Cultivation and transfection conditions of experiment 1 (Ambr® 15).

Process parameters	Set points
Stirring speed (rpm)	600,800
pH	6.9; 7.1; 7.3
Constant parameters	30% DO; 15 mL fill volume; 37 °C cultivation temperature
Transfection conditions	1 µg DNA + 2 µl PEIpro®/10 ⁶ cells; plasmid ratio: 5:2.5:1.5:1 (GOI:gag-pol:VSV-G:rev)

These two factors were evaluated as to their contribution to the lentivirus titer yield (Figure 2). According to Figure 2A, cells generally grew better in the Ambr® 15 than in the control shake flask culture. However, large differences in the fold expansion can also be observed between the different conditions tested in the Ambr® 15. The viability of the cells cultured at pH 7.3 drops significantly at day 2 (not shown) which explains the reduced lentivirus titer at this pH. Overall, transient transfection with PEIpro® did not impact cell viability (data not shown). Generally, it was observed that higher pH values and stirring speeds yielded improved cell growth (Figure 2A), but these factors negatively correlate to LV particle titer as seen in the DoE model (Figure 2B). This may imply that a 2-step approach could be beneficial with a shift of the main process parameters when transitioning from the growth phase to the production phase at the time of the plasmid transfection.

We also identified that the lentivirus titer is higher in the Ambr® 15 vessels than in the positive control shake flask. According to Figure 2A we could clearly identify optimal culture conditions for lentivirus production in the Ambr® 15 microbioreactor. A stirring speed of 600 rpm and a pH between 6.9 and 7.1 yielded the highest lentivirus titer ($8.8 \times 10^9 - 9.8 \times 10^9$ VP/mL). This trend was confirmed with the DoE model. According to the response contour plot a clear trend of an increasing lentiviral particle titer could be observed with a decreasing stirring speed and a peak lentiviral titer was observed at pH culture values between 7.0 and 7.1. Furthermore, it can be concluded that the culture pH

and the stirring speed are critical process parameters that have a significant effect on the viral vector production.

OPTIMIZATION OF TRANSFECTION WITH PEIPRO® FOR LENTIVIRUS PRODUCTION IN AMBR® 15

To optimize the transient transfection process, we used PEIpro® transfection reagent in the Ambr® 15 microbioreactor system. PEIpro® benefits from extensive research development that make this unique PEI-based transfection reagent optimal for lentivirus production in adherent and suspension systems system [13]. This is in part due to its unique ability to efficiently condense several plasmid DNA and deliver them into HEK-293 cells for production of full lentivirus particles. We optimized four parameters that could have an impact on the success of transient transfection and therefore on virus production: the viral production medium, DNA amount, ratio of DNA/transfection reagent and ratio of plasmids.

Through a DoE approach we were able to screen all these parameters in one cultivation run, thanks to the screening capabilities of the Ambr® 15 microbioreactor system [5-7]. Recapitulated in Table 2 are the ranges tested for each parameter, based on previous experience and manufacturers protocols. A D-optimal design with triplicate center points was chosen, leading to 23 different conditions / vessels.

According to Figure 3A, the LV particle titer is higher in some of the Ambr® 15

vessels than in the shake flask positive control. Depending on the conditions used for transfection, extreme differences in titers between the culture vessels can be observed, indicating that the factors analyzed indeed have a significant effect on lentivirus productivity. At optimal conditions, a viral titer of 2.1×10^{11} VP/mL and a specific productivity, meaning the particle titer per cells at the time of transfection, of 1.3×10^5 VP/cell was obtained which was higher than the titer obtained in the reference shake flask (1.4×10^{11} VP/mL and 5.9×10^4 VP/cell).

The infectious viral titer of the best condition in the Ambr® 15 seems to be equal to the positive control shake flask, however,


when comparing the specific productivity of the lentivirus, meaning the infectious titer per cells at the time of transfection, this value is significantly higher for the Ambr® 15 vessel (20.13 for the Ambr® 15 vs. 13.66 TU/cell for the reference shake flask).

Furthermore, when comparing the highest viral particle titer obtained in this lentivirus production run with the highest one from the first experiment, we see another twentyfold increase in viral particles through this optimization.

After obtaining the LV particle titer for all culture vessels, we analyzed the DoE model of this screening experiment with the MODDE® software. Our results lead to a good

► **FIGURE 2**

(A) Lentivirus titer and HEK-293T/17 SF cells fold expansion during the optimization of the culture conditions in the Ambr® 15. (B) Results graphs from the analysis of the DoE model with MODDE®.

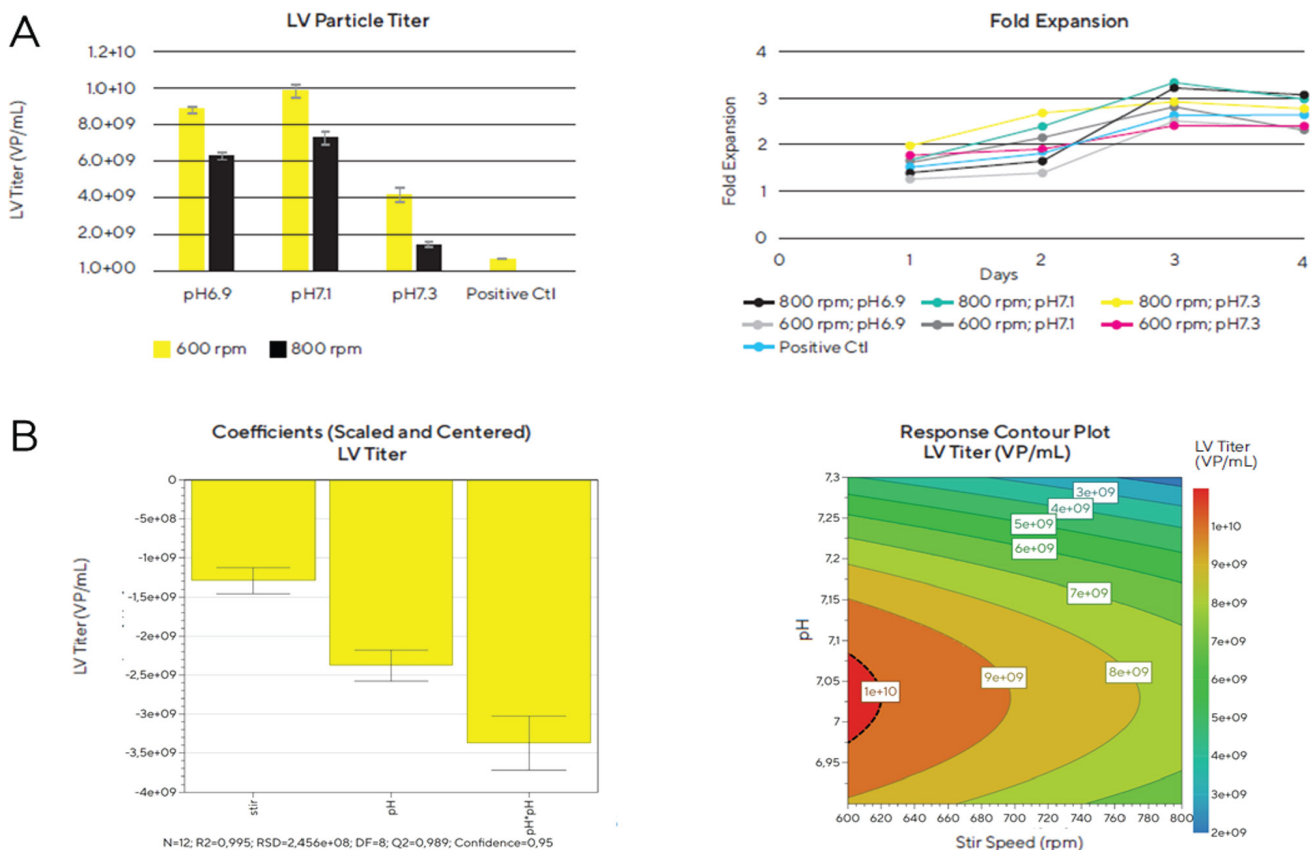

(A) Shown are mean values of duplicate vessels with standard deviation. Positive control = standard shake flask culture. (B) All conditions have been analyzed in duplicates. The experimental design (full factorial design with two replicates of each condition) was created with MODDE® software. Model coefficient factors on the left (stir speed and pH) and their impact on the process readout lentivirus particle titer. In the right, a response contour plot with LV particle titer profile.

▶ TABLE 2
Cultivation and transfection conditions of experiment 2 in Ambr® 15.

Process parameters	Set points/ranges
Constant parameters	600 rpm; pH limits 6.9–7.1; 30% DO; 15 mL fill volume; 37 °C cultivation temperature
Viral production medium	Freestyle293; SFM4Transfx-293
DNA amount	0.5 – 4µg DNA/10 ⁶ cells (at transfection)
Ratio PEIpro®:DNA	1:1 – 4:1
Plasmid ratio: GOI:gag-pol:VSV-G:rev	5:2.5:1:1; 5:1:2.5:1; 5:1:1:2.5

modeling of the DoE for the optimization process. Furthermore, when we plotted the results in a response contour plot, we could clearly see an optimal spot of the lentivirus particle titer when the ratio of PEIpro® to DNA was high and the DNA amount per 10⁶ cells was low. We were able to identify several factors that have a significant effect on the lentivirus titer during the transfection process. For example, the ratio of the amount of PEIpro® to DNA, the usage of Freestyle 293 medium and the plasmid ratio of 5:2.5:1:1 positively correlate to LV particle titer. However, the amount of DNA, the usage of SFM4Transfx 293 medium and a plasmid ratio of 5:1:2.5:1 negatively correlate with the lentivirus titer. The plasmid encoding the gene of interest (GOI) was always present in excess due to the larger size of this plasmid with the GOI to improve virus particle production capacity. Reducing the VSV-G amounts while increasing gag-pol amounts provided a good balance between proteins needed for virus replication and proteins involved in the LV enveloped particles formation.

Therefore, with only one production run we were able to identify optimal set points of key factors which led to a twentyfold increase in lentivirus titer compared to the initial protocol used in experiment 1.

According to the optimizer function of the DoE software MODDE® our new optimal conditions for the transfection step during lentivirus production are:

DNA amount: 0.5 µg/10⁶ cells; ratio PEIpro®:DNA: 4:1; viral production medium: Freestyle 293; plasmid ratio: 5:2.5:1:1 (GOI:gag-pol:VSV-G:rev)

UPSCALING LENTIVIRUS PRODUCTION PROCESS IN AMBR® 250 MODULAR

After having optimized the transient transfection process step and cultivation parameters for lentivirus production, we tested the feasibility of upscaling the optimized process to the Ambr® 250 Modular. We furthermore aimed to optimize the gas flow rate and stir direction which is enabled in this bioreactor system with automated processing. These parameters could potentially have a significant impact in the viral titer due to its sensitivity to externally applied forces (i.e., shear forces) and the dependence of the viral titer on the cell viability.


The culture parameter set points and transfection conditions are listed in Table 3. Due to the limited scalability of the Ambr® 15 system, the optimal stirring speed for the production process in the Ambr® 250 Modular was identified by running a separate experiment which included testing of four different stir speeds in the Ambr® 250 Modular.

According to the results shown in Figure 4, the lentiviral particle titer was significantly higher in vessel/ condition 3 of the bioreactor system which corresponds to a maximum gas flow rate of 0.5 mL/min and a down stirring direction of the impellers. In this vessel, a viral titer of 2.1 x 10¹¹ VP/mL and a specific productivity of 1.3 x 10⁵ VP/cell was obtained which was higher than the titer obtained in the reference shake flask (8.9 x 10¹⁰ VP/mL and 3.7 x 10⁴ VP/cell).

The highest viral titer was obtained with the lowest gas flow rate and we could prove that the gas flow rate and the stir direction have a significant effect on lentiviral particle

► **FIGURE 3**

(A) Lentiviral particle and infectious titer obtained in a screening experiment to optimize the transfection process with the Ambr® 15. (B) Analysis of the DoE model of the transfection conditions screening experiment.

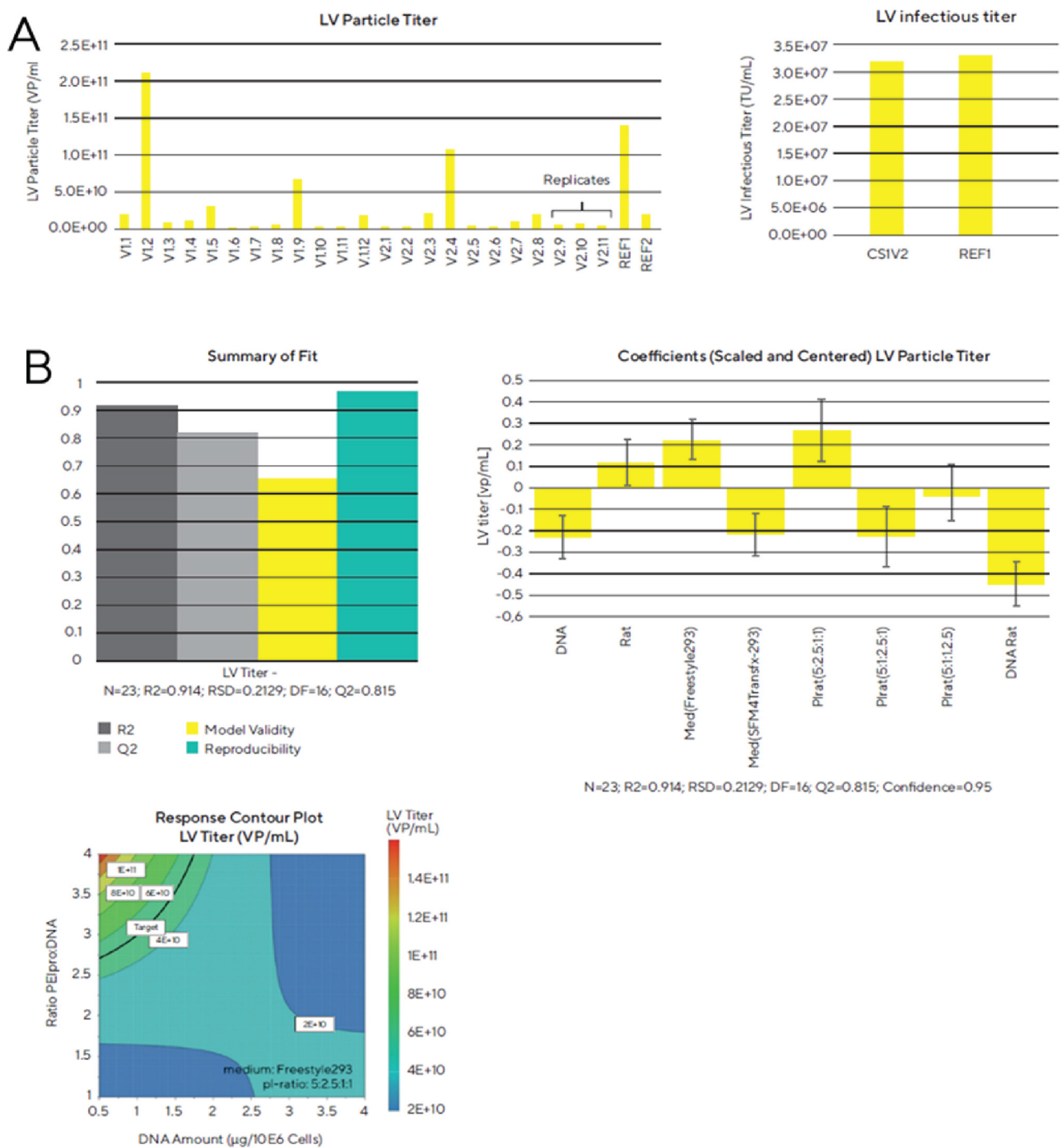

(REF1/2=shake flask control with each of the tested media; each bioreactor vessel represents a different transfection condition as defined by the DoE layout)

titer, indicating a negative impact of high gas flow rates on lentiviral titer. Varying flow rates and stir directions also led to differences

in fold expansion of the cells during lentivirus production. Even though the viral particle titer was very high in vessel 3, we also observed

TABLE 3
Gas flow rates and stir directions of experiment 3 (Ambr® 250 Modular).

Process parameters	Set points	
Constant parameters	400 rpm; pH 7.1; 30% DO; 250 mL fill volume; 37 °C cultivation temperature	
Transfection conditions	0.5 µg DNA + 2µl PEIpro®/10 ⁶ cells; plasmid ratio: 5:2.5:1:1 (GOI:gag-pol:VSV-G:rev)	
Vessel/condition	Gas flow rates (air/mix) [mL/min]	Stir direction
1	0.1–0.5	Up
2	0.1–2.5	Up
3	0.1–0.5	Down
4	0.1–2.5	Down
All:	CO ₂ and O ₂ added flow: 0–5	

a good growth rate of the cells in this bioreactor. In general, the cells cultured in vessel 1 and 3 showed a better growth profile than vessel 2 and 4, indicating that the gas flow rate has a major impact on cell growth and viability. In this experiment we were able to show that the optimized lentiviral vector production protocol is scalable to larger bioreactor volumes, and that the gas flow rate had a significant impact on lentiviral titer. With the most optimal condition, we were able to obtain a lentiviral particle titer of 2.1×10^{11} VP/mL and a specific productivity of 1.3×10^5 VP/cell in Ambr® 250 Modular which was consistent with the optimized LV production in Ambr® 15 (2.1×10^{11} VP/mL and 1.3×10^5 VP/cell). These results confirm that production of lentiviral vectors can directly

be scalable from the Ambr® 15 microbioreactor system to a larger stirred bioreactor system without loss in yield.


CONCLUSION

This study demonstrates that the Ambr® 15 microbioreactor system in combination with the DoE software MODDE® enables a systematic investigation of critical process parameters and rapid, high throughput process optimization in a reduced time. Optimization of cultivation parameters and of the transfection process is critical to substantially improve lentiviral titer.

The results prove that the transition from shake flask to a scalable stirred bioreactor

FIGURE 4

Lentiviral particle titer and fold expansions of the HEK-293T/17 SF cells obtained in an upscaling experiment to optimize the gas flow rate and stir direction with the Ambr® 250 Modular.

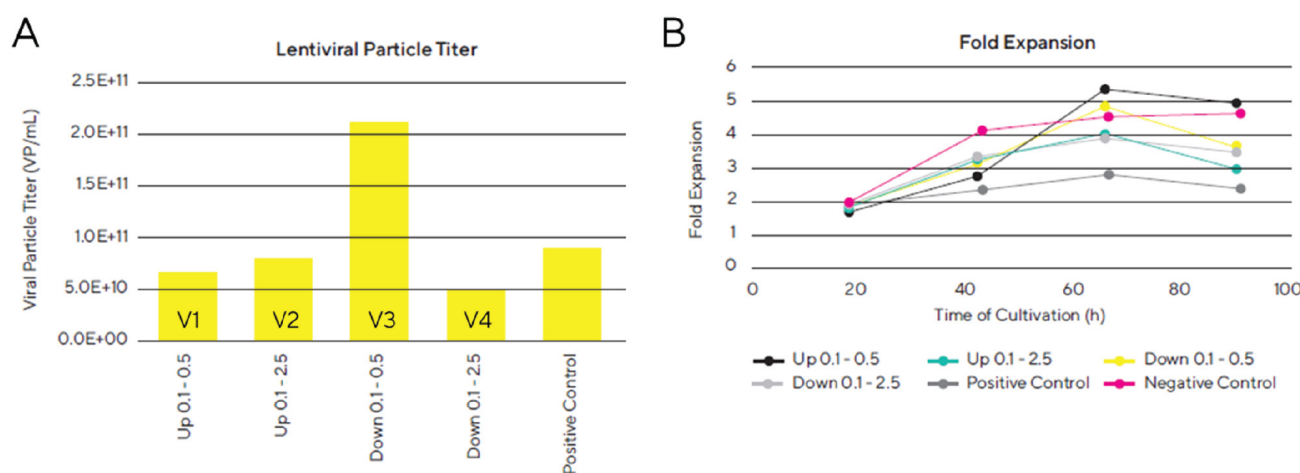

system can be facilitated and lead to further improved lentiviral titers. Optimization can directly be ensured with the right set of tools: the automated Ambr® 15 microbioreactor system combined with a scalable transfection process with PEIpro® transfection reagent. Scale-up of the process is simplified by relying on the Ambr® 250 Modular. The Ambr® 15 correlated well with the Ambr® 250 Modular results and provides good basis for further scale-up to larger stirred bioreactors as shown previously for a mAb process [11].

The outcome of such a study is designed to help manufacturers gain important process knowledge on the parameters that need to be controlled to set up a robust and predictable lentivirus production process that supports scaling to much larger scales (up to 2k L) for GMP manufacturing at commercial scale.

METHODS

Lentivirus production

Third generation lentivirus was produced by transient transfection of suspension HEK-293T/17 SF cells (ATCC #ACS-4500) in a stirred bioreactor (either Ambr® 15, Ambr® 250 Modular).

The cells had been passaged at least twice before starting the lentivirus production.

Freestyle 293 media (Thermo Fisher Scientific) was filled into the respective bioreactors on day 0 and process parameter control was initiated. Later, on day 0, the cells were seeded into the bioreactor at a final VCD of 1×10^6 cells/mL. A lentivirus production in a 125 mL shake flask was prepared equally to the bioreactor (positive control).

After 24 h, transfection of a CD19-CAR encoding transfer plasmid and three lentiviral helper plasmids (Aldevron) are performed using PEIpro® DNA transfection reagent (Polypius Transfection). A defined amount of DNA (sum of all four plasmids) per 1×10^6 cells is diluted in Freestyle 293 medium at a certain

plasmid ratio (the volume is 1:20 of the final culture volume). In a separate reaction tube a defined volume of PEIpro® per 1×10^6 cells is diluted in Freestyle 293 medium (the volume is 1:20 of the final culture volume). Diluted PEIpro® is added to the diluted DNA, gently mixed and incubated for 15 min at room temperature. The mixture is added dropwise to the cells. A negative control without using a transfection reagent is prepared and treated equally (cells are cultured in a 125 mL shake flask).

On the next day, i.e. 18 h after transfection, anti-clumping reagent (1:500 (v/v), Thermo Fisher Scientific) and 10 mM sodium butyrate (Sigma) are added.

LV was harvested 72 h post transfection. Before harvesting, the virus suspension was treated with 10 U/mL DENARASE® (c-Lectta) for 1 h for digestion of nucleic acids.

Lentivirus quality control & analysis

As a primary readout on virus concentration, we performed a p24-ELISA, that measures lentivirus-associated p24 protein, to determine the total viral particle titer. The assay was performed according to the manufacturer's protocol (Cell Biolabs). The assay's accuracy was determined to be below 8 % CV.

Due to the nature of typical infectious titer assays, being very laborious and giving low sample throughput, we decided to primarily run a particle titration assay (p24-ELISA) and only determine the infectious titer of selected samples based on the viral particle titer results. Still, viral particle titers allow for observation of overall effects of factors on the lentivirus production process.

A flow cytometry-based assay was performed to determine the infectious lentiviral titer by transducing adherent HEK-293T cells with the lentiviral supernatants [12,14].

The HEK-293T/17 SF cell density and viability were measured with a Cedex HiRes instrument (Roche).

REFERENCES

1. Anguela XM, High KA. Entering the Modern Era of Gene Therapy. *Ann. Rev. Med.* 2019; 70(1): 273–88.
2. Levine BL, Miskin J, Wonnacott K, Keir C. Global Manufacturing of CAR T Cell Therapy. *Mol. Ther. Methods Clin. Dev.* 2017; 31; 4: 92–101.
3. Young H, Philip Probert P. Development and scale-up of suspension culture processes for viral vector manufacturing: challenges and considerations. *Cell Gene Ther. Ins.* 2020; 6(1): 149–57.
4. Warnock JN, Merten OW, Al-Rubeai M. Cell culture processes for the production of viral vectors for gene therapy purposes. *Cytotechnology* 2006; 50(1–3): 141–62.
5. Krakowiak J, Liu Q, Nascimento-Brooks L. Rapid development of viral vector production processes. *BioProcess Int.* 2021; 19(4).
6. Moses S, Manahan M, Ambrogelly A, Ling WLW. Assessment of AMBR as a model for high-throughput cell culture process development strategy. *Adv. Biosci. Biotechnol.* 2012; 3: 918–27.
7. Shahid Rameez, Sigma S. Mostafa, Christopher Miller, Shukla AA. High-throughput miniaturized bioreactors for cell culture process development: Reproducibility, scalability, and control. *Biotechnol. Prog.* 2014; 30(3): 718–27.
8. Manahan M, Nelson M, Cacciatore JJ, Weng J, Xu S, Pollard J. Scale-down model qualification of ambr® 250 high-throughput mini-bioreactor system for two commercial-scale mAb processes. *Biotechnol. Prog.* 2019; 35(6): e2870.
9. Janakiraman V, Kwiatkowski C, Kshirsagar R, Ryll T, Huang YM. Application of high-throughput mini-bioreactor system for systematic scale-down modeling, process characterization, and control strategy development. *Biotechnol. Prog.* 2015; 31(6): 1623–32.
10. Sandner V, Pybus LP, McCreath G, Glassey J. Scale-Down Model Development in ambr systems: An Industrial Perspective. *Biotechnol. J.* 2019; 14(4): e1700766.
11. Ruhl S, de Almeida N, Carpio M, Rupperecht J, Greller G, Matuszczyk JC. A Rapid, Low-Risk Approach Process Transfer of Biologics from Development to Manufacturing Scale. *BioProcess Int.* 2020; 18(5).
12. Labisch JJ, Bollmann F, Wolff MW, Pflanz K. A new simplified clarification approach for lentiviral vectors using diatomaceous earth improves throughput and safe handling. *J. Biotechnol.* 2020; 326: 11–20.
13. Nyamay'antu A, Kedinger V, Erbacher P. Addressing scaling-up limitations: optimized PEI-mediated production of clinical grade viral vectors. *Cell Gene Ther. Ins.* 2018; 4(2): 71–9.
14. Kutner RH, Zhang XY, Jakob Reiser J. Production, concentration and titration of pseudotyped HIV-1-based lentiviral vectors. *Nat. Protoc.* 2009; 4(4): 495–505.

AFFILIATIONS

Diana Riethmüller

Sartorius Stedim Biotech GmbH,
Göttingen, Germany

Dr Alengo Nyamay'antu

Polyplus-transfection, Illkirch, France

Dr Franziska Bollmann

Author for correspondence:
Sartorius Stedim Biotech GmbH,
Göttingen, Germany
franziska.bollmann@sartorius.com



AUTHORSHIP & CONFLICT OF INTEREST

Contributions: All named authors take responsibility for the integrity of the work as a whole, and have given their approval for this version to be published.

Acknowledgements: The authors acknowledge Polyplus Transfections for support in publishing the study and Lara Nascimento-Brooks, Market Entry Strategy Manager at Sartorius, for support with the manuscript and its publication.

Disclosure and potential conflicts of interest: *The authors declare that they have no conflicts of interest.*

Funding declaration: *The authors received no financial support for the research, authorship and/or publication of this article.*


ARTICLE & COPYRIGHT INFORMATION

Copyright: Published by Cell and Gene Therapy Insights under Creative Commons License Deed CC BY NC ND 4.0 which allows anyone to copy, distribute, and transmit the article provided it is properly attributed in the manner specified below. No commercial use without permission.

Attribution: Copyright © 2021 *Name INTL*. Published by Cell and Gene Therapy Insights under Creative Commons License Deed CC BY NC ND 4.0.

Article source: Invited; externally peer reviewed.

Submitted for peer review: May 12 2021; **Revised manuscript received:** Jun 18 2021; **Publication date:** *date*.

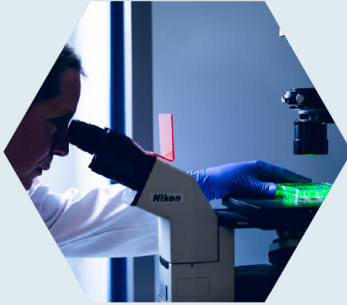

First to provide GMP grade reagent for viral vector manufacturing



Supplier of FectoVIR®-AAV:
Next generation of transfection reagent for recombinant AAV virus production in suspension cells

With 4 transfection product lines, Polyplus-transfection is the leading biotechnology company that supports Gene and Cell therapy, biologics manufacturing and life science research with innovative nucleic acid transfection solutions.



Viral vectors for Gene & Cell Therapy



Life science research



Protein and mAb production



Non viral *in vivo* transfection

20
years

20 years expertise in transfection solutions



Available worldwide

www.polyplus-transfection.com

Contact us: